Analyzing T-RFLP Files Using Peak Scanner

Background:

Once upon a time, and a very good time it was, there was a simple way to analyze T-RFLP data. First one used GeneScan to generate the peak files, then one followed with a couple of wonderful scripts written by Dr. Zaid Abdo at the University of Idaho which prepared the data for downstream clustering analysis. However, as time went on, GeneScan got older and ABI ceased to support it. This would still have been fine had ABI also made sure that chromatogram files generated on their newer machines could not be read by GeneScan. Let me emphasize that last point again since I've wasted an obscene amount of time trying to search for the backward compatibility that should be there: Chromatogram files generated on any ABI sequencer that is less that 4 or 5 years cannot be read by GeneScan. It is worth noting; however, that there is a conversion program that was made by ABI, but they will not give it out. Of course even if you had the conversion program GeneScan does not play nicely with XP. Fortunately, ABI has just released a free program called Peak Scanner which fills in many of the gaps left by not supporting GeneScan.

What you will need:

First you will need to go the ABI website and download Peak Scanner (type in Peak Scanner into the search if needed). Peak Scanner is a Windows only program (known to work under XP SP2, I don't know about Vista). If you are using an Intel Mac Peak Scanner runs beautifully in Parallels.

Peak Scanner Setup:

Once you have downloaded Peak Scanner you have got to set it up to be able to analyze the T-RFLP data.

1) Start Peak Scanner and you will be presented with the interface pictured in Figure 1.

	📖 Mordoor – Parallels Desktop		
New Open Save Close Edit. Print Export Ar	chive Analysis	👔 🗸 Help	Preferences
To begin, select from the following options:			Ċ
New Project Select this option to start a new project, which will require you to add samples, a analysis method and size standard to each sample, and then analyze the sample and sizing information.	ossign an to view data Settings\All Users\App	Open Project ts from the table below or use the 'Browse' function to select a pr the projects listed in the table below are located at "C:\Documents ication Data\Applied Biosystems\Peak Scanner\Application Data".	oject and
	Project Name	Type Last Modified	
Start New Project	Open Project:	Open Selected Project	
	http://www.appliedbiosystems.com		
i Peak Scanner		6) 9:19 PM
Click inside OS Window to capture mouse		H D 0 0	9 0 0 18 //

Figure 1

2) Next click on the New button in the upper left hand corner and select New Size Standard. You should see a window pop up as depicted in Figure 2. Name the

k Scanner					. 2 🛛	
e Standard Editor						
					U v 🕾	
Ste Standard Name:	20X1000				nep meetroos	
Site standard rand.						e.,
Description:	R0X-1000					
Dye Color:	Red 💌					
					of function to celect a moject	(FR
Enter sizes in the field	below separated by a comma se	pace, or return then click the	"Add Site(s)ss" bullon to add them to th	e current size standard definition	cated at "C:\Documents and	
		,			nner'(Application Data".	
Enter new Size Standa	rd definition: (e.g. 11.0, 34.2, 55	5)	Current Size Standard definition:	Delete Selected Sizes	1	
			50.0	^		-
			75.0			
			100.0			
			125.0	=		
			300.0			
			250.0			
		Add Size(s) >>	300.0			
			350.0			
			400.0			
			450.0			
			475.0			
			500.0	~		
		Save As	Save	Cancel		
					browse	
	Start New Project			Open Selected Project		
		bite	Www.emlachionstam.com			
		000				

size standard and set the fluorophor color. Then enter in the sizes (for ROX-1000 it should be 50, 75, 100, 125, 150, 200, 250, 300, 350, 400, 450, 475, 500, 550, 600, 650, 700, 750, 800, 850, 900, 950, 1000). Click on Save or Save As.



3) Select New again only this time select New Analysis Method. You should see a window open up like in Figure 3. Name the method (T-RFLP Analysis or something to that effect). Next

	💭 Mordoor – Parallels Desktop		click on Peak
sis Method Editor Name: T-RFLP Analysis Method Description: Pask smoothing Range Baselining Minimum peak heights	Torwsef function set Scamer (Apple) Modified	Preferences	Smoothing, Range, Peak Characteristics, Size Calling Method, Baselin ing, and Minimu Peak Heights ar fill in the values shown in Figures 4-9 respectively.
Revert to Defaults Revert	Save As Save Cancel	Figure 4	Figure 3
	Heavy The default is set to "None". If the peaks appear jagged, applying smoothing will refine the displayed peaks.		

Peak smoothing Range		Name	Blue	Green	Yellow	Red	Orange	
Peak characteristics	1	Minimum Peak Half Width	1	1	1	1	1	
Size calling method Baselining	2	Polynomial Degree	3	3	3	3	3	
Minimum peak heights	3	Peak Window Size	15	15	15	15	15	
	4	Peak Slope Start	0.0	0.0	0.0	0.0	0.0	
	5	Peak Slope Stop	0.0	0.0	0.0	0.0	0.0	
	These settings determine peak characteristics. Minimum Peak Half Width: any peak narrower than this setting (in scans) will not be called as a peak. Polynomial: changing this number will increase/decrease sensitivity of peak morphology,							
	Th Mir pe Po res	ese settings determine peak nimum Peak Half Width: any p ak. ynomial: changing this numb pectively.	characte beak nar er will ine	eristics. rower than t crease/decre	his setting (in so ase sensitivity c	ans) will no	t be called as phology,	a

Figure 6

Peak Detection Quality Flags Peak smoothing Range Peak characteristics Stee colling method Baselining Minimum peak heights	 Global Southern Least Square Local Southern Cubic Spline 	Figure 7
	Select a size calling method. Local Southern is the recommended method.	

Peak Detection	Quality Flags			
Peak smoothin Range Peak characte Size calling me	ig ristics thod	Bartha Uta dan Gara - Fi		
Baselining Minimum peak	heights	baseline window Size: 51		Figure 8
		Baseline Window Size determines the size of the window to be used for baselining. Increasing the window size will decrease the baselining effect.		
			~	
		<u>x</u>	2	

Peak smoothing Range Peak characteristics Size calling method	Baseline Window Size: 51		
Baselining Minimum peak heights			Figure 9
	Baseline Window Size determines the size of the window to be used for baselining. Increasing the window size will decrease the baselining effect.		
		~	

Next click on the Quality Flags tab and select Sizing Quality and fill in the values as shown in Figure 10.

ing quality		
	Sizing Quality	
	Pass Range: From 0.75 to 1.0	
	Low Quality Range: From 0.0 to 0.25	Figure 1
	Assume Emelancy From (ep.), 30.0 TO (ep.); 000.0	
	These settings will determine the pass/fail status of the size standard.	
	If the sizing quality fails in the Low Quality Range, no peaks will be sized.	
	The range specified in 'Assume Linearity' indicates that the size standard fragments falling within the specified range, migrate linearly.	
		~
		and the second se

And Save the method. Now you are ready to analyze some data.

4) Click on the Start New Project button and you will be presented with a Project Window as in Figure 11.



Click on the Add Files button and navigate to where the chromatogram files are located. If you are using Parallels and the files were downloaded in the Mac environment don't worry, just drag the folder that contains the files to the Windows desktop and the folder will be copied over.

Figure 11

Now select the Chromatogram files you want to analyze, click on the Add Selected Files button then click on OK. The project window should now look something like what is

🔹 - 🧀 - 🔚 - 🤤 ew Open Save Cl	📕 📝 🗸 🍋 🗸 📸	l → rt	Archive	- 🕨 - Analysis					🕜 🗸	et al la constant de
Project > Untitled	Display: 16 Plot	_								
	To Add Country and an the									
Setup	To Add Samples, click on the	s woo r	nes button	below or urag and un	op itsa nies into c	ane application v	window.			
Add Samples To Project	🗊 Samples View	Add File	Remove	Files Size Standard	🖂 Analysis Me	sthod 🖂	Show/Hide Samples:	Show All	× 🗔 🛙	10
Set Applyois Developmentary	Sample Explorer <<	1	Status	Sample Name	Sample Type	Size Stan	dard Analysis M	ethod SQI	Offscale	Quality
Set Analysis Parameters	IIII Untitled	1	A.	38908_A01	sample	~	*	~		
(CONTRACTOR OF CONTRACTOR OF C	😐 🔛 3486-Chromatogram	r 2	2	38908_A02	sample	~	~	~		
analyze		3	л?	38908_A03	sample	~	~	~		
		4	1.º	38908_A04	sample	~	N	~		
Quality Control		5	1. L	38908_A05	sample	~	~	~		
Edit Size Standards		6	2	38908_A06	sample	~	~	~		
		7	H.	38908_A07	sample	~	~	~		
		8	22	38908_A08	sample	~	~	~		
Review Data		9	10	38908_A09	sample	~	×	M		
View		10	25	38908_A10	sample	~	M	<u> </u>		
I abal Daske		11	25	38908_A11	sample	M	×	<u> </u>		
		<					_			>
Edit Peaks		碧	Not Ready	for Analysis 🛛 🦉 Rea	ady for Analysis	Analyzed	File Corrupted	File Locked	File Not Fou	und 🗸
2					(2	24) problem(s) found.			
Resources		Sam	ple File Nar	ne		Description			Pro	blem Id
Now Look ale Method		3486	5.A01.fsa			Size standa	ard not specified		3	
New Analysis Method		3486	5.A01.fsa			Analysis me	ethod not specified		7	
New Size Standard		3486	AUS.Isa			Size standa	ard not specified		3	
Manage Analysis Methods		3488	A10 fee			Sine shands	ethod not specified		2	
		3400	A10.fsa			Dize standa	and not specified		3	
		3484	AD4 fca			Size stands	and not specified		3	
Applied		3400	new made			Jace Stallue	and not specified		5	

shown in Figure 12. Now don't worry about the error messages they will be fixed shortly.

Go to the Size Standard column of the first row and click on the tab. Select the Rox-1000 that you created earlier (or other desired size standard). Now click on the Size Standard column so that the whole column is selected and type Ctrl-d.

Figure 12

This should fill down the column with the same selection. If it doesn't work deselect the column, reselect it and try again. This happens sometimes and is part of this being a free (and a bit buggy) program. Next click on the tab in the first row of the Analysis Method column and select the T-RFLP method that you created. Select the whole column and fill down as before. Now everything should be set and no more errors should be reported. Finally, click on the Analyze Files.

5) Getting your peak files. You should see a window that looks something like what is shown in Figure 13.



From here on in the order in which you do things can make a huge difference. Also which subwindow you have selected when you go to do a function makes a big difference so be aware of where you are clicking. So in order do the follow:

1) in the graph subwindow click on the rainbow square so that all of the colors are deselected. Next click on the blue square so that the blue channel is selected (Figure 14)

Not Ready for Analy	sis 🦉 Ready for Ana Data: Ani	ilysis 🍌 Analyzed 🔋	File Corrupted	ile Locked 👔 File Nol 💥 🔟 Mode:	t Found 🛛 Pass Basepairs 👽 🛄	Fail Check	-
] 3486.A01.fsa 0 200	400	600 800	1000 1	.200 1400	>>	Analyzed Data Dye Legend	
400		-*****	***		-12000	1.0 3486.A01.fsa	
200 -					-4000 0 <		Figure
			Show A	All Peaks	Label Selecte	d Peaks 📄 👩 🖻	-
Sizing Table View			SHOW: SHOTT				

2) Select all of the files from the file list subwindow (Ctrl-a)

3) From the graph subwindow click on the Edit Plot Settings button

This will bring up a new window although for some reason it almost always brings it up behind the main window. So look to the task bar and you should see that one of them is titled Plot Settings, click on that to bring the settings window to the front. In the settings window click on the Labels Tab and then click on the Label Peaks button as shown in Figure 15.



4) Next click on the Edit Table Settings button in the top left of the bottom subwindow as shown in Figure 16.

Sizing	Table View			Show: Show All Peaks	~	Label Selected Peaks		-
	Dye Color	Dye/Sample Peak	Sample File Name	Size	Height	Area in Point	Area in I	^
1		B, 1	3486.A01.fsa	50.3395	3	16	1	
2		B, 2	3486.A01.fsa	51.6123	6	49	4	
3		В, З	3486.A01.fsa	52.8844	6	28	2	
4		B, 4	3486.A01.fsa	53.9864	7	66	5	
5		B, 5	3486.A01.fsa	55.1726	6	32	2	
1				54 0500				-

Figure 16

This will bring up the Table Settings window. Make the Columns Selected on the right hand side of the window look like Figure 17. Click on Apply to apply this formatting. Before closing the window click on Save As to save this formatting. In the future you will just be able to select this method from the drop down menu and then hit Apply to format the table. Keep in mind that in order for downstream data analysis to work it is vitally important that this formatting be exactly as you see it.

lect the columns to be displayed. Adjust Available Columns to Display	the order displayed in the table u	sing i	<i>the move up/move a</i> Columns	<i>lown controls.</i> Selected	
Dye Color		[Column Name	Sort Order	
oampie File Name Δrea in RP		2	Begin Point		
Begin BP		3	Size	~	
End Point		4	Height	~	Figure ⁻
End BP	Add Selected >>	5	Area in Point	~	
Width in Point		6	Data Point	~	
Width in BP					
User Comments					
User Edit	<< Remove Selected				
Project Name	<< Remove All				
Run Name					
Run Date					
Run Duration					

6) In the bottom subwindow click on the Show drop down menu and select Show Labeled Peaks (Figure 18). This will confine the data in the tables to just the blue channel.

	Table View			Show:	Show Labeled Peaks	V Lab	el Selected Peaks 🛛 🔚 🕜 🖵	
	Dye/Sample Peak	Begin Point	Size	Height	Show All Peaks Show Selected Peaks	pata Point	^	
1	B, 1	1795	50.3395	3	Show Labeled Peaks	1801		
2	B, 2	1806	51.6123	6	49	1816		
3	В, З	1823	52.8844	6	28	1831		
4	B, 4	1836	53.9864	7	66	1844		
5	B, 5	1852	55.1726	6	32	1858		
6	B, 6	1864	56.3583	4	30	1872		

Set up for file export is now complete.

7) Select the first file in the file list. Now click on the Sizing Table View tab of the bottom subwindow. Now click on the Export button at the top of the window and select Export. Save the files in a new folder. You should name the files sequentially and something simple like A1 or M1. Repeat this step with every file. Remember the order to this because if you forget to click on the Sizing Table View tab you will export the wrong thing. When you have exported all of the files just quit Peak Scanner. Don't bother saving the project because yet another bug of the program is that you can never get a saved project to work properly. Remember it's free though.

8) Now drag the file that has all of the Peak Files onto the Mac Desktop or wherever for downstream analysis.

Troubleshooting:

I didn't write the program, but I've used this enough to field some questions. So if you have questions about this tutorial or basic Peak Scanner problems that are not covered here, then I can be reached at jre@umich.edu.